


## APPLICATION NOTE

No. 8

### Nexygen DF setup

KRD 9-9-20

#### Configuring

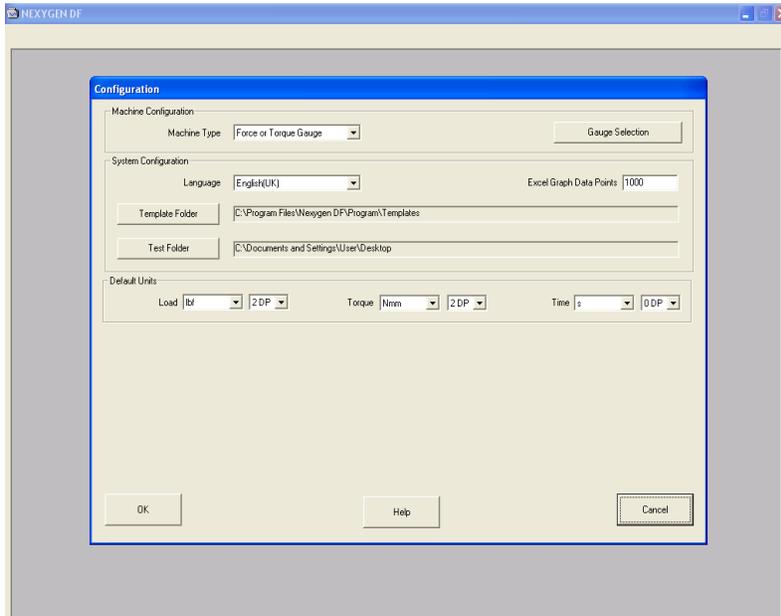

Select gage type and properties.

Change Excel Data Points

CAUTION: Do not change Template or Test folder location.


If Template Folder is changed to the desk top, no other location may be specified in the program until:

- Nexygen DF in uninstalled completely and then reinstalled

If Test Folder location is changed to desktop follow the above procedure OR

- Temporarily change Template Folder to `C:/`, then when pressing Test Folder enter is quickly pressed.

#### Creating a Test

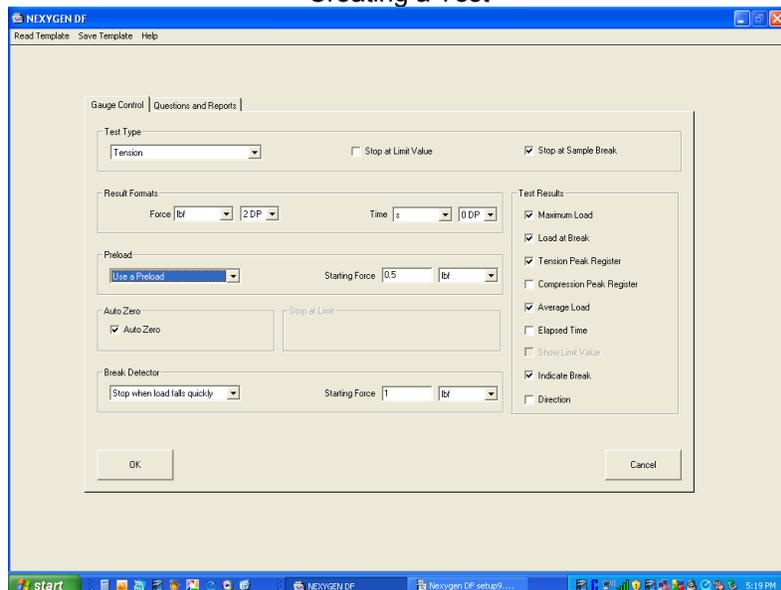

Set Test Type

Set how test will end. Selecting Stop At Break will enable Break parameters. This portion will stop the graphing automatically

Set Preload if desired. This portion will allow the graph to automatically start graphing

Select the results desired.

## APPLICATION NOTE

### Setting up Export, Questions and Reports

Add in any questions you want to have included in you test.

Enter the Report Setup Locations

Enter EXCEL Export location.


If you will be using this setup profile again, save it as a template.

Note: No changes to this page can be made once a test has been saved

When you click ok, this screen will be displayed.

CAUTION: File name must be at least 4 characters long. If it is shorter, the program will lock up later in use.

## APPLICATION NOTE

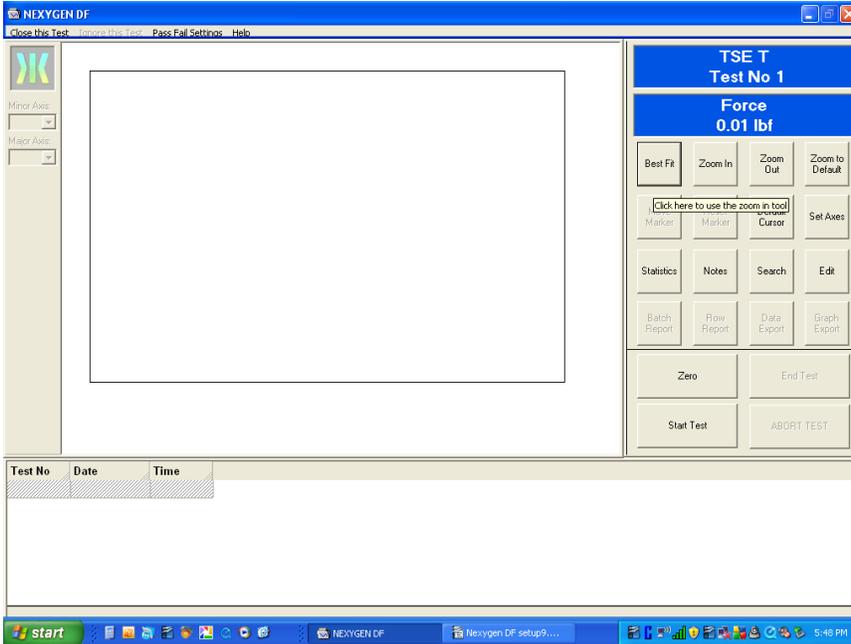


After saving file, you are ready to test.

Press Start Test to start graphing. If Preload was selected in the setup, graphing will not start until the threshold force is achieved.

To stop the test without collecting data, press Abort Test.

If no End Test parameters were set up, press End Test to stop Graphing

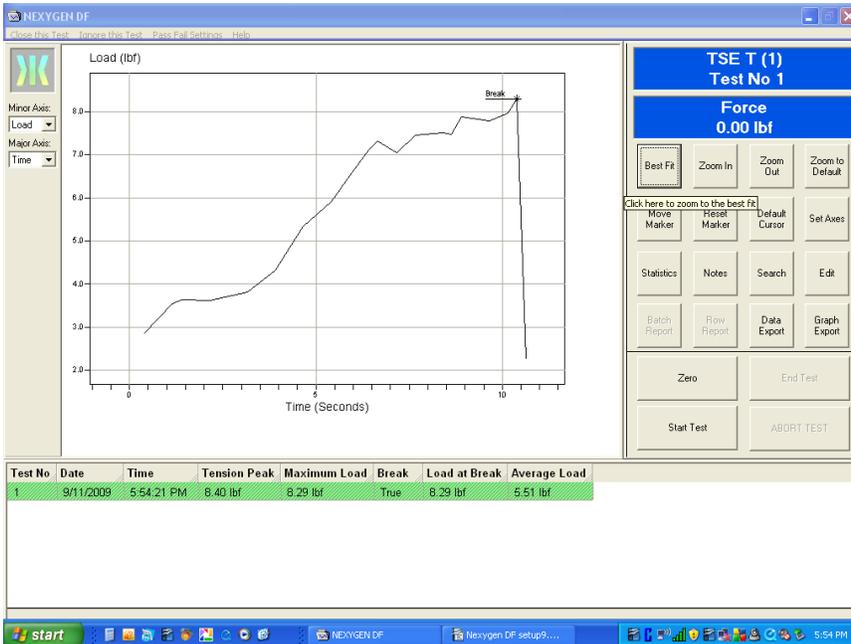

After the test has ended data or graph may be exported by pressing the Data Export or Graph Export button.

Note: these buttons will not be highlighted if an Excel file was not setup earlier.

Caution: it is not possible to setup data or graph export after the test is run.